

GSE

Economy Magnetic Stirrer

User Manual

What's in the box:

- Magnetic Stirrer
- Power Adapter
- User Manual

PRODUCT DESCRIPTION

The GSE Magnetic Stirrer from Globe Scientific is an excellent choice for most basic stirring applications. The unit comes with integrated retort stand to hold electrochemistry probes, temperature probes and more. A spill-resistant chassis and universal voltage adapters are included.


SAFETY PRECAUTIONS

The detailed precautions described in this User Manual were carefully developed in an attempt to cover all the possible risks. Please read carefully before operating this stirrer.

- Ensure that only trained staff works with the stirrer.
- Never use the stirrer in any manner not specified in these instructions.
- If any abnormal condition occurs during operation, please stop use immediately and contact Globe Scientific.
- **Caution – Magnetism!** Effects of the magnetic field must be considered (e.g. data storage media, cardiac pacemakers, etc.).

OPERATION

Place the GSE stirrer on a level, firm surface and confirm the power supplied matches the power supply's rated voltage. When ready to operate, rotate the speed adjustment knob to control the GSE stirring speed, clockwise increases the speed, counterclockwise decreases the speed. The GSE stirrer can be adjusted from 300rpm to 2000rpm.

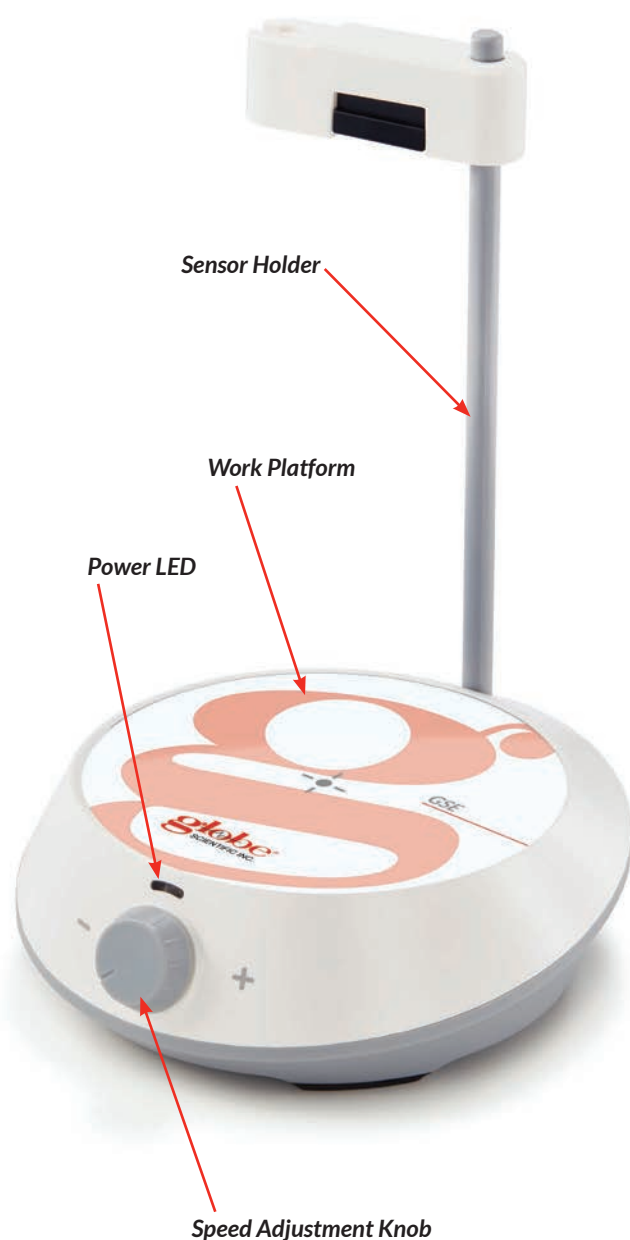

Continued ...

GSE

Economy Magnetic Stirrer

User Manual

GSE Specifications	
Voltage	100-240 VAC
Frequency	50/60 Hz
Power	1.2W
Adjustment	Scale
Speed Range	300 – 2000rpm
Max. Volume (H ₂ O)	1.5L
Max. Stir Bar Size	35mm
Stir Platform	120mm dia.
Stir Platform Material	PET plastic
Dimension (WxDxH)	145 x 150 x 50mm (5.7 x 5.9 x 1.9in)
Weight	0.3kg (0.66lbs)
Operating Temperature Range	5-40°C (41-104°F)
Permissible Humidity	80% non-condensing
Protection Class	IP21
Warranty	2 Years

WARRANTY

This stirrer is guaranteed for two years from the date of delivery provided it has been operated and maintained. Spills, corrosion and misuse are not covered and will void this warranty.

SERVICE:

In the unlikely event you require assistance of any kind, please contact Globe Scientific immediately.



www.globescientific.com



Customer Service, Sales and Technical Support:

Phone: 1 (800) 394-4562 • 1 (201) 599-1400 ext. 2

Fax: 1 (201) 599-1406 • E-mail: support@globescientific.com