


GTR-2-AS / Adjustable Speed Tube Rocker Quick Start Guide

What's in the box:

- Quick Start Guide
- Tube Rocker
- Power Cables (1 each: US plug, UK plug, EU plug)
- 10 Position Silicone Tray Inserts

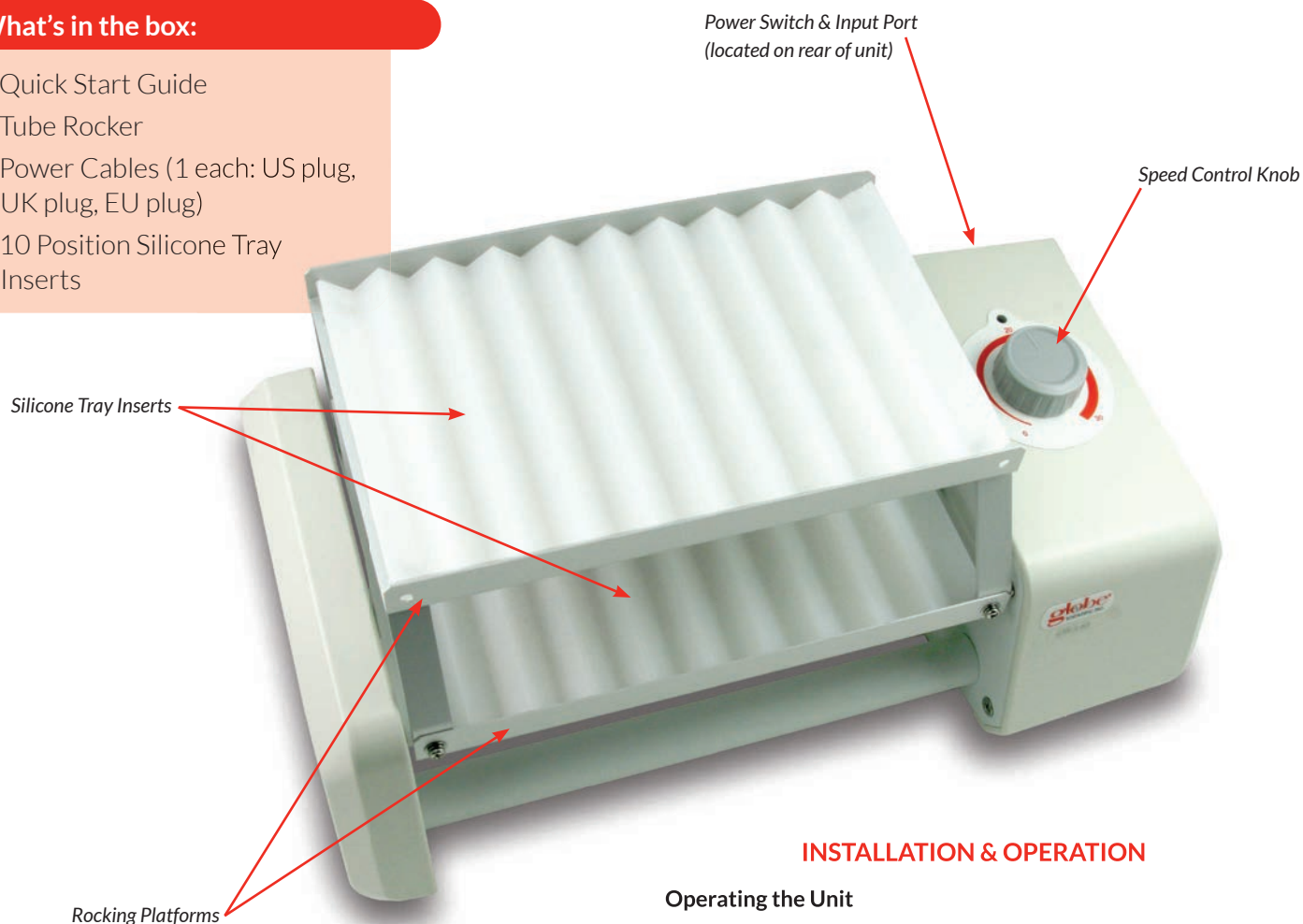

IMPORTANT: The tube rocker must not be overloaded or damage may occur. Please do not exceed 1kg (2.2 lbs).

INSTALLATION & OPERATION

Operating the Unit

- Carefully unpack and place the unit on a stable non-slip surface.
- Allow approximately 4 inches surrounding the unit for safe operation.
- Connect the power adaptor using the correct plug for your power supply.
- Turn power switch on.
- Rotate the speed control knob to the desired speed setting.



Customer Service, Sales and Technical Support:

Phone: 1 (800) 394-4562 • 1 (201) 599-1400
Fax: 1 (201) 599-1406 • E-mail: mail@globescientific.com