

GTR-2-FS Tube Rocker Quick Start Guide

What's in the box:

- Tube Rocker
- Power Cable(s)
- Quick Start Guide

OPERATION

- Simply switch the main power to the rocker ON or OFF. When the unit is turned ON, it will rock at a fixed speed of 25rpm.

Power Switch



IMPORTANT:

- The tube rocker must not be overloaded or damage may occur. Please do not load the unit in excess of 7lbs (3.2kg).

www.globescientific.com

globe[®]
SCIENTIFIC INC.

Customer Service, Sales and Technical Support:

Phone: 1 (800) 394-4562 • 1 (201) 599-1400

Fax: 1 (201) 599-1406 • E-mail: mail@globescientific.com