

GTR-AVS / Variable Speed Analog Tube Roller Quick Start Guide

What's in the box:

- Quick Start Guide
- Tube Roller
- Power Cables (3 each)
- Extra Fuse



CONTROL PANEL



Power Indicator

LED light will glow green when unit is powered on.

Speed Control Knob

Rotate clockwise to start operation and increase speed. Rotate counterclockwise to decrease speed and stop operation .

On / Off Switch

Toggles power to the unit on and off.

IMPORTANT: The tube roller must not be overloaded or damage may occur. Do not load the unit in excess of 8.8lbs (4kg).

NOTE: When handling the tube roller, always lift the unit by the front and rear housing only, never by the connecting rods.

For the full User Manual, please visit www.globescientific.com



Customer Service, Sales and Technical Support:

Phone: 1 (800) 394-4562 • 1 (201) 599-1400

Fax: 1 (201) 599-1406 • E-mail: mail@globescientific.com