

GTR-DVS / Variable Speed Digital Tube Roller Quick Start Guide

What's in the box:

- Quick Start Guide
- Tube Roller
- Power Cables (3 each)
- Extra Fuse



CONTROL PANEL

Speed / Timer Indicator

Green LED indicates the LCD screen currently displays speed. Red LED indicates the LCD screen currently displays time.

LCD Display

Displays the unit status and all set values.

On / Off Switch

Toggles power to the unit on and off.



Power / Overload Indicator

Green LED indicates power on. Steady red LED indicates overload protection is enabled. Flashing red LED indicates overload and roller operation is stopped.

Mode Switch Button

Press to switch between speed and time.

Speed / Timer Control Knob

Rotate to set speed and time. Push to start/stop.

IMPORTANT: The tube roller must not be overloaded or damage may occur. Do not load the unit in excess of 8.8lbs (4kg).

NOTE: When handling the tube roller, always lift the unit by the front and rear housing only, never by the connecting rods.

For the full User Manual, please visit www.globescientific.com



Customer Service, Sales and Technical Support:

Phone: 1 (800) 394-4562 • 1 (201) 599-1400

Fax: 1 (201) 599-1406 • E-mail: mail@globescientific.com