

GVM-FS

Vortex Mixer

Quick Start Guide

What's in the box:

- Vortex Mixer
- Quick Start Guide

CONTINUOUS OPERATION

- Place the vortex mixer on a stable and safe surface and plug in the power cord.
- Press the toggle switch **left** to ON to enable Continuous mode. The Power LED will light.
- Push the toggle switch right lightly to the middle position to turn unit OFF when mixing is complete.

TOUCH OPERATION

- Place the vortex mixer on a stable and safe surface and plug in the power cord.
- Press the toggle switch **right** to Touch to enable Touch mode.
- When a test tube is pressed into the mixing head vertically, the vortex mixer will begin to mix and the Power LED will light.
- Push the toggle switch left lightly to the middle position to turn unit OFF.



On / Off Switch

Toggles power between continuous mode, touch mode and off.



NOTE:

- Place vessels and accessories as vertical as possible into the center of mixing head for proper mixing.
- Do not run continuous operation mode indefinitely.

Continued ...


GVM-FS

Vortex Mixer

Quick Start Guide



CHANGING THE ATTACHMENT

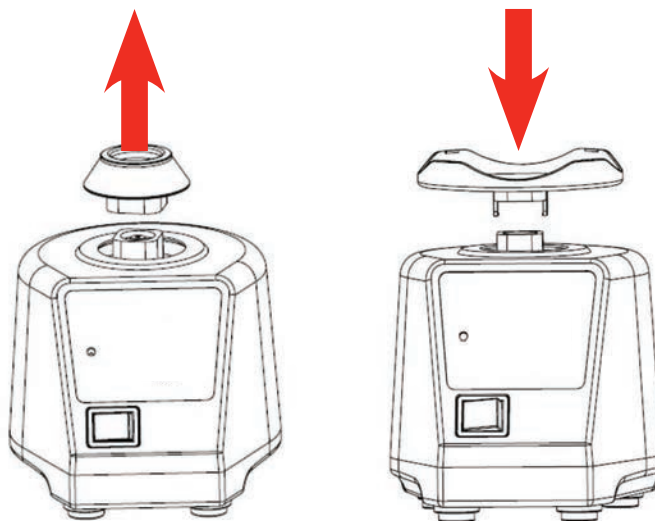

ACCESSORY COMPATIBILITY CHART

| ATTACHMENT | SUITABLE FOR: | |
|-----------------------------------|---------------|-----------------|
| | Touch Mode | Continuous Mode |
| GVM-AS-CUP | YES | YES |
| GVM-AS-ROD and GVM-AS-ADAPT48 | YES | NO |
| GVM-AS-ROD and GVM-AS-ADAPT18 | YES | NO |
| GVM-AS-ROD and GVM-AS-ADAPT12 | YES | NO |
| GVM-AS-ROD and GVM-AS-ADAPT8-16 | YES | NO |
| GVM-AS-ROD and GVM-AS-ADAPT8-20 | YES | NO |
| GVM-AS-PLATE and GVM-AS-PAD | YES | NO |
| GVM-AS-PLATE and GVM-AS-ADAPT48 | NO | YES |
| GVM-AS-PLATE and GVM-AS-ADAPT18 | NO | YES |
| GVM-AS-PLATE and GVM-AS-ADAPT12 | NO | YES |
| GVM-AS-PLATE and GVM-AS-ADAPT8-16 | NO | YES |
| GVM-AS-PLATE and GVM-AS-ADAPT8-20 | NO | YES |

www.globescientific.com



Customer Service, Sales and Technical Support:

Phone: 1 (800) 394-4562 • 1 (201) 599-1400

Fax: 1 (201) 599-1406 • E-mail: mail@globescientific.com