

GVM-MINI

Vortex Mixer

Quick Start Guide

What's in the box:

- Vortex Mixer
- Power Adaptor
- Quick Start Guide

TOUCH OPERATION

- Place the vortex mixer on a stable and safe surface and plug in the power cord.
- Plug the power adaptor in the appropriate voltage wall outlet as indicated on the adaptor.
- Ensure that the label indicated the correct voltage before connecting the device to power supply.
- Switch power ON and observe whether the power indicator (LED) is lit.
- The test tube should be pressed into the center of the mixing head with as much force as necessary to allow the mixer to work steadily.
- Release the force applied on the mixing head to stop the mixer.

NOTE:

- Place vessels and accessories as vertical as possible into the center of mixing head for proper mixing.
- Do not run continuously for an indefinite amount of time.



www.globescientific.com

globe[®]
SCIENTIFIC INC.

Customer Service, Sales and Technical Support:

Phone: 1 (800) 394-4562 • 1 (201) 599-1400

Fax: 1 (201) 599-1406 • E-mail: mail@globescientific.com